

SEPABEADS™ SP700

SEPABEADS™ SP700 is highly porous styrenic adsorbents. It has highest surface area in SEPABEADS™ series and a narrower pore size distribution than DIAION™ HP20. It can be used for various application.

SEPABEADS™ SP700 is characterized by:

- >> Higher surface area and unique pore size distribution
- >> High chemical and physical stability
- >> Excellent pressure/flow characteristics
- >> Excellent batch-to-batch reproducibly
- >> Wide application

Physical and chemical properties

Grade Name	DIAION™ SP700	
Bead form	Spherical, porous	
Matrix	Poly divinylbenzene / ethylvinylbenzene	
Chemical Structure	$ \begin{array}{c} \text{---CH}_2\text{---CH---CH}_2\text{---CH---} \\ \qquad \qquad \\ \text{C}_6\text{H}_4 \qquad \text{C}_6\text{H}_4 \\ \qquad \qquad \\ \text{---CH---CH}_2\text{---} \qquad \text{CH}_2\text{CH}_3 \end{array} $	
Whole Bead Count	-	95 min.
Shipping Density*	g/L	690
Water content	%	60 - 70
Particle Size Distribution thr. 250 μm	%	5 max.
Effective size	mm	0.25 min.
Uniformity Coefficient	-	1.6 max.
Particle Density*	g/mL	1.02
Specific Surface Area	m ² /g	1100 min.
Pore Volume*	mL/g	2.2
Pore Radius*	Å	90

Note : properties with a mark "*" are referential data.

Swelling ratio in various solvents

Methanol	1.04
Ethanol	1.05
2-Propanol	1.07
Acetone	1.07
Toluene	1.05
Acetonitrile	1.05
Water	1.00



Pore size distribution

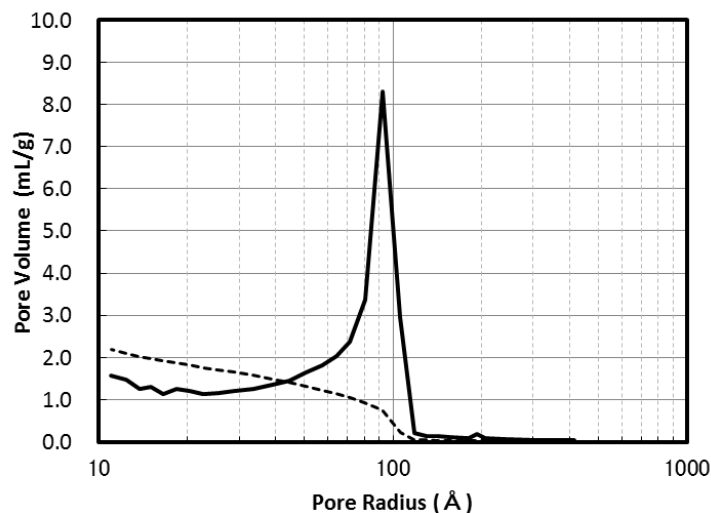


Fig. 1 Pore size distribution of SP700

Recommended Operating Conditions

Maximum Operating Temperature	°C	130
Operating pH Range		0 - 14
Minimum Bed Depth	mm	800
Flow rate	BV/h	Loading 0.5 - 5
	BV/h	Displacement 0.5 - 2
	BV/h	Regeneration 0.5 - 2
	BV/h	Rince 1 - 5
Regenerant		
	Organic solvents for hydrophobic compounds	
	Bases for acidic compounds	
	Acids for basic compounds	
	Buffer solution for pH sensitive compounds	
	Water for an ionic solution	
	Hot steam for volatile compounds	



Hydraulic Characteristics

The approximate pressure drop at various temperatures and flow rates for each meter of bed depth of SEPABEADS™ SP700 resin in normal down flow operation is shown in the graph below.

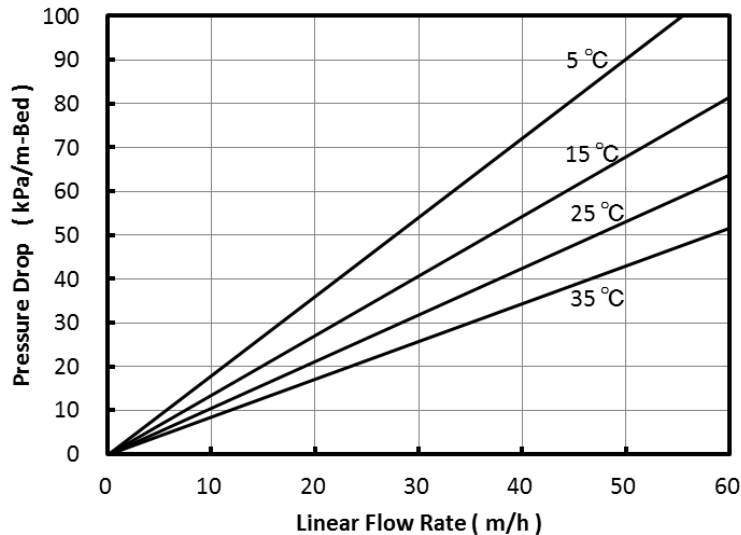


Fig. 2 Pressure Drop of SP700

FDA status

SEPABEADS™ SP700 may be used to process food and beverage products and isolate specialized food additives as intended and such used may be said to fully comply with the Federal Food, Drug, and Cosmetic Act.

Applications

- Purification of juices
- Removal of naringin and other bittering agents
- Purification of small peptides, oligonucleotides and proteins
- Adsorption of vitamins, antibiotics, enzymes, steroids and other substance from fermentation solutions
- Decolorization and purification of various chemicals

Notice

This information are given in good faith but without warranty, and this also applies where proprietary rights of third parties are involved. The application, use and processing of our products are beyond our control and therefore your own responsibility.

