



RAPID EPS Instruction Manual

Thank you for purchasing our RAPID EPS.
Please read this manual carefully before use and save this manual for your future use.

Please note that product specifications and packings may change without prior notice.

Announcement for your use with ease.

【Purpose of use】

This product can be used for research purposes only.

【Notes for use】

Handle this product carefully when unpacking from the bag.

Do not process this product in autoclave treatment.


【Storage and disposal of product】

Store at room temperature and in a place not exposed to direct sunlight. To dispose the product after use, comply with local laws and ordinances, and process properly according to the regulations of your facility.

Instructions

*Please wear gloves to handle the product.

1

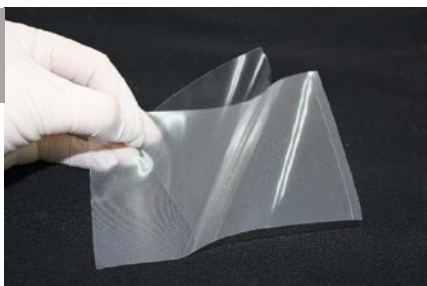

Remove the release liner from the slit along the edge and expose the adhesive surface.

2



Adhere the adhesive surface on a well plate. After pasting, press the surface with hands or a spatula, etc. and **adhere it firmly.**



1040 1st Avenue
Suite # 330
New York, NY 10022
Phone: 212-204-0075
Email: info@pyvot.tech
Web: www.pyvot.tech

Catalog Code	IUREPS001 PV03-1003
Product Name	RAPID EPS
Bonding method	Pressure sensitive adhesive
Materials	polyolefin
Operating temperature limit	-80°C~80°C
Size	W 80 mm × L 122 mm
Quantity	100sheets/Box