

SUPER SlitSeal

Instruction Manual

Thank you for purchasing our SUPER SlitSeal.

Please read this manual carefully before use and save this manual for your future use.

Please note that product specifications and packings may change without prior notice.

Announcement for your use with ease.

【Purpose of use】

This product can be used for research purposes only.

【Notes for use】

Handle this product carefully when unpacking from the bag.


Do not process this product in autoclave treatment.

【Storage and disposal of product】


Store at room temperature and in a place not exposed to direct sunlight. To dispose the product after use, comply with local laws and ordinances, and process properly according to the regulations of your facility

Instructions

*Please wear gloves before use.

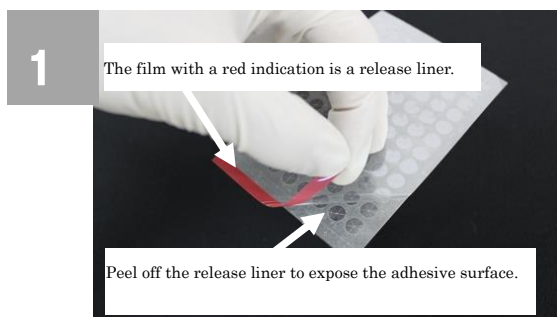

Remove the release liner with a red indication and expose the adhesive surface.

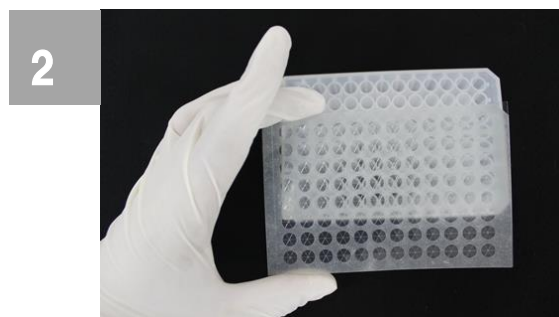

Paste the SUPER SlitSeal on the 96-well plate and make sure holes of SUPER SlitSeal are correctly positioned on each well spot.



Remove the protective film before sampling or dispensing. If used without removing the protective film, the needle or the tip can not puncture the film. If the protective film stays on and is left at 100°C, the edge of the protective film may be curved.



1040 1st Avenue
Suite # 330
New York, NY 10022
Phone: 212-204-0075
Email: info@pyvot.tech
Web: www.pyvot.tech

Catalog Code	IURSS-S96-801- 22 PV03-1002
Product name	SUPER Slit Seal
Bonding method	Pressure sensitive adhesive
Materials	PET, Silicone
Operating temperature limit	-80°C~100°C *
Size	W 80 mm × L 122 mm
Quantity	100 sheets/Box

* The upper limit of the operating temperature is specified as 100 ° C, and this is clarified by the evaluation test carried at 100 ° C for 30 minutes.